

Программное обеспечение SET

Технические характеристики

По вопросам продаж и поддержки обращайтесь:

Алматы (7273)495-231	Калининград (4012)72-03-81	Омск (3812)21-46-40	Сыктывкар (8212)25-95-17
Ангарск (3955)60-70-56	Калуга (4842)92-23-67	Орел (4862)44-53-42	Тамбов (4752)50-40-97
Архангельск (8182)63-90-72	Кемерово (3842)65-04-62	Оренбург (3532)37-68-04	Тверь (4822)63-31-35
Астрахань (8512)99-46-04	Киров (8332)68-02-04	Пенза (8412)22-31-16	Тольятти (8482)63-91-07
Барнаул (3852)73-04-60	Коломна (4966)23-41-49	Петрозаводск (8142)55-98-37	Томск (3822)98-41-53
Белгород (4722)40-23-64	Кострома (4942)77-07-48	Псков (8112)59-10-37	Тула (4872)33-79-87
Благовещенск (4162)22-76-07	Краснодар (861)203-40-90	Пермь (342)205-81-47	Тюмень (3452)66-21-18
Брянск (4832)59-03-52	Красноярск (391)204-63-61	Ростов-на-Дону (863)308-18-15	Ульяновск (8422)24-23-59
Владивосток (423)249-28-31	Курск (4712)77-13-04	Рязань (4912)46-61-64	Улан-Удэ (3012)59-97-51
Владикавказ (8672)28-90-48	Курган (3522)50-90-47	Самара (846)206-03-16	Уфа (347)229-48-12
Владимир (4922)49-43-18	Липецк (4742)52-20-81	Саранск (8342)22-96-24	Хабаровск (4212)92-98-04
Волгоград (844)278-03-48	Магнитогорск (3519)55-03-13	Санкт-Петербург (812)309-46-40	Чебоксары (8352)28-53-07
Вологда (8172)26-41-59	Москва (495)268-04-70	Саратов (845)249-38-78	Челябинск (351)202-03-61
Воронеж (473)204-51-73	Мурманск (8152)59-64-93	Севастополь (8692)22-31-93	Череповец (8202)49-02-64
Екатеринбург (343)384-55-89	Набережные Челны (8552)20-53-41	Симферополь (3652)67-13-56	Чита (3022)38-34-83
Иваново (4932)77-34-06	Нижний Новгород (831)429-08-12	Смоленск (4812)29-41-54	Якутск (4112)23-90-97
Ижевск (3412)26-03-58	Новокузнецк (3843)20-46-81	Сочи (862)225-72-31	Ярославль (4852)69-52-93
Иркутск (395)279-98-46	Ноябрьск (3496)41-32-12	Ставрополь (8652)20-65-13	
Казань (843)206-01-48	Новосибирск (383)227-86-73	Сургут (3462)77-98-35	
Россия +7(495)268-04-70	Киргизия +996(312)-96-26-47	Казахстан +7(7172)727-132	

KERN SET-01

ET OS à KERN EasyTouch operating system with basic functions ET Lab-Set + ET Industry-Set
(weighing + laboratory package)

Software EasyTouch

OS Operating System SET-01

Category

Brand: Accessories KERN

Category: Software

Product Group: Software

Packaging & shipping

Delivery: 24 h

Shipping method: Parcel service

Piktogramme

STANDARD



FACTORY



KERN SET-04

ET Density à Density determination function

Software EasyTouch

Density SET-04

Category

Brand: Accessories KERN

Category: Software

Product Group: Software

Packaging & shipping

Delivery: 24 h

Shipping method: Parcel service

Piktogramme

STANDARD



OPTION



KERN SET-05

ET Dynamic â Animal and dynamic weighing function (e.g. animal weighing mode)

Software EasyTouch

Dynamic SET-05

Category

Brand: Accessories KERN

Category: Software

Product Group: Software

Packaging & shipping

Delivery: 24 h

Shipping method: Parcel service

Piktogramme

STANDARD



OPTION



KERN SET-06

ET Classify à Classification function

Software EasyTouch

Classify SET-06

Category

Brand: Accessories KERN

Category: Software

Product Group: Software

Packaging & shipping

Delivery: 24 h

Shipping method: Parcel service

Piktogramme

STANDARD



OPTION



KERN SET-07

ET Variable à Free Variable Function

Software EasyTouch

Variable SET-07

Category

Brand: Accessories KERN

Category: Software

Product Group: Software

Packaging & shipping

Delivery: 24 h

Shipping method: Parcel service

Piktogramme

STANDARD




OPTION



KERN SET-08

ET Take-out â The take-out function enables negativeweighing out of a source container

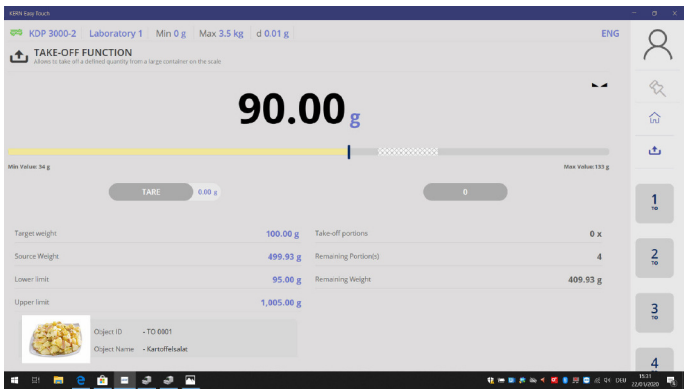

Category	
Brand:	Accessories KERN
Category:	Software
Product Group:	Software
Packaging & shipping	
Delivery:	24 h
Shipping method:	Parcel service
Piktogramme	

STANDARD



OPTION



KERN SET-10

ET Save-Server à Weight storage and Server storage function

Software EasyTouch

Save-Server SET-10

Category

Brand: Accessories KERN

Category: Software

Product Group: Software

Packaging & shipping

Delivery: 24 h

Shipping method: Parcel service

Piktogramme

STANDARD



OPTION



KERN SET-11

ET Safety à Precision Safety function

Software EasyTouch

Safety SET-11

Category

Brand: Accessories KERN

Category: Software

Product Group: Software

Packaging & shipping

Delivery: 24 h

Shipping method: Parcel service

Piktogramme

STANDARD



OPTION



KERN SET-14

ET Print â User-customised print format function

Piece Counting

Master Object ID	C00012
Master Object Name	Fixing Element #17 649
Dynamic Object id	FEX #17e9578
Dynamic Object name	Fixing Element Sample
Quantity	1429 pcs
Net Result	200.10 g
Tare weight	3.57 g
Gross Weight	203.67 g
Reference weight	0.140 g
Date	2020-01-22
Time	17:30:00
User	admin

Used Device	KDP 3000-2
Serial Number	W1805832
Last Adjustment	2020-01-22
Internal Code	Laboratory 1

KERN & Sohn GmbH, Ziegelei 1, 72336, Balingen, Germany
Tel: 0049 7433 9933, E-mail: info@kern-sohn.com

Category

Brand:	KERN
Category:	Software
Product Group:	Software

Packaging & shipping

Delivery:	24 h
Gross weight:	0 mg
Net weight:	0 mg
Shipping method:	Parcel service

Piktogramme

STANDARD




OPTION



KERN SET-16

ET Batch and Statistics â Batch measuring function with statistical analysis and batch print function

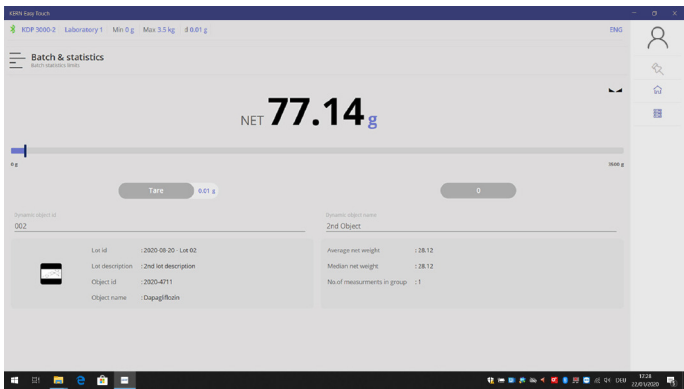

Category

Brand:	KERN
Category:	Software
Product Group:	Software

Packaging & shipping

Delivery:	24 h
Shipping method:	Parcel service

Piktogramme

STANDARD




OPTION



KERN SET-21

ET Formulation à Recipe function

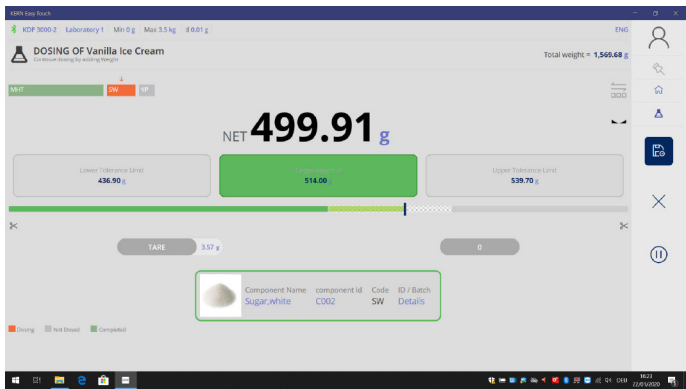

Category

Brand:	Accessories KERN
Category:	Software
Product Group:	Software

Packaging & shipping

Delivery:	24 h
Shipping method:	Parcel service

Piktogramme

STANDARD




OPTION



KERN SET-22

ET Difference à Difference weighing e.g. the growth of cell cultures

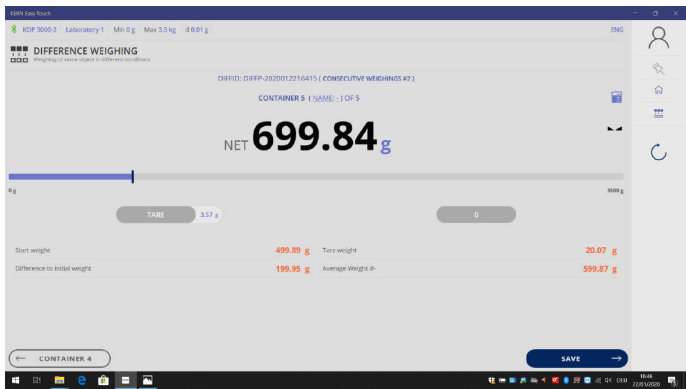

Category	
Brand:	Accessories KERN
Category:	Software
Product Group:	Software
Packaging & shipping	
Delivery:	24 h
Shipping method:	Parcel service
Piktogramme	

STANDARD



OPTION



Software SET-31

ET Tolerance – Tolerance weighing function

NEW

Software EasyTouch **Tolerance** **SET-31**

- Note: the required basis is SET-01 ET OS (basic package). SET-31 is part of SET-03 Premium Industrial package with “Piece-counting function” and “Target-counting function” at a bargain price
- The tolerance-weighing App allows rapid, simple as well as complex checkweighing procedures. For this, the user specifies a lower and an upper weight limit. Using the EasyTouch system, a range of objects can be weighed and the system indicates whether that particular weight is within the defined limits or is outside those limits. This function in many applications, including sorting, portion division as well as order picking
- You can enter the limits very quickly in EasyTouch Tolerance using the touchscreen or by entering them in the appropriate field on the PC. ^[1] The limits can be entered in two ways: either just with the lower and upper limit (absolute limit) ^[3] or using a target value with an upper and lower permitted deviation ^[2]. The permitted deviations can be entered either in “g” (as an absolute value) or in “%” of the target value (as a relative value) ^[5]
- ^[6] Central master data memory: Checkweighing objects can be stored in the memory of the system with a target weight, as well as lower and upper limits. In this way these limits do not have to constantly be entered again, but can be easily recalled from the memory. In the master data memory you can also store a possible tare value for the typical packaging, box or container which is typically used for the object and which will then be deducted automatically from the weighing result

- **4** ID security: It offers the possibility of storing each weighed and stored weighing result with a unique ID number (Dynamic Object ID) and an ID name (Dynamic Object Name). Storage can be triggered either semi-automatically or fully automatically and this will always happen after the weight has been removed from the balance and then a load has been applied. This means that the user does not have to press any keys to save the weights and can work efficiently
- Colour bar graph: Through the colour bar graph, the user is shown quickly and clearly, whether the weighing result is below, within or above the tolerance range. The red “I” result marking also shows the user where exactly the weight is within the tolerance range.
- Full mode: In full mode, the complete weighing range of the screen is shown in colour, with the relevant colour for the tolerance weighing result. In this way the user has an even clearer indication of the relevant tolerance result
- **9** Batch ID: In addition for “ID security” of each individual object, a “Batch ID” can be assigned for a full test batch. This batch ID number is saved with each individual memory. In this way, you guarantee that later all stored results in the dynamic data memory can be identified with this batch ID
- **10** Tolerance function in pieces: It is possible to operate the checkweighing function in pieces instead of g/kg. To do this, enter the individual weight of the objects to be counted (reference weight) as well as the target quantity and the upper and lower limit quantities
- Tolerance function in percent: It is possible to operate the checkweighing function as a percentage instead of g/kg. To do this, enter the target value as well as the upper and lower values in percent.
- Watch the product film

Technical data

- Licensing: One license can be operated on up to four terminal devices (PC, laptop, tablet) at the same time, working independently
- User: You can store as many users as you need in one license
- Balances: You can store and operate as many balances as you need in one license
- Communication balances-terminal device: The balance(s) can communicate using serial connections, USB, Bluetooth, Ethernet or WiFi with the PC, laptop or tablet

STANDARD



OPTION




ALL MODELS

Model

SET-31

KERN SET-32

ET Count à Piece-counting function

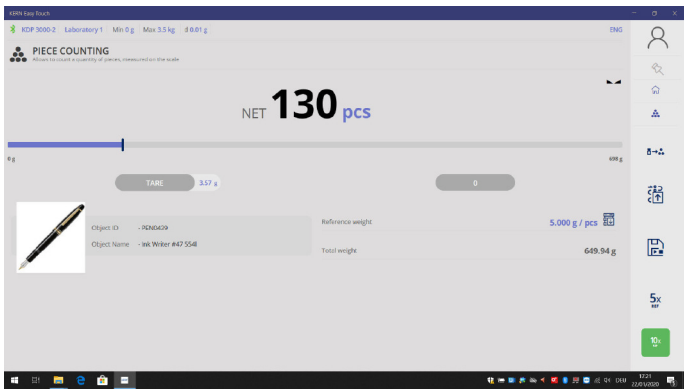

Category

Brand:	Accessories KERN
Category:	Software
Product Group:	Software

Packaging & shipping

Delivery:	24 h
Shipping method:	Parcel service

Piktogramme

STANDARD




OPTION



KERN SET-33

ET Target-count â Target-counting function

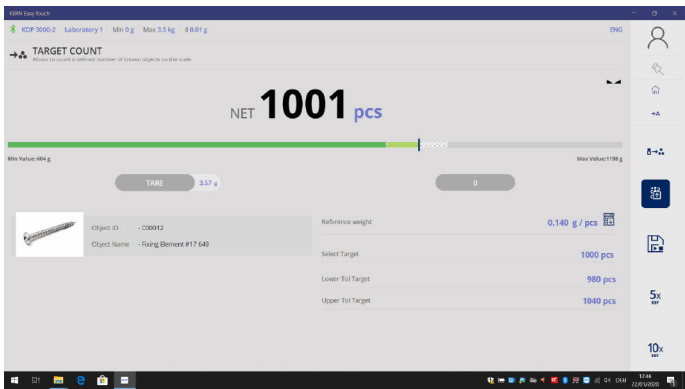

Category

Brand:	Accessories KERN
Category:	Software
Product Group:	Software

Packaging & shipping

Delivery:	24 h
Shipping method:	Parcel service

Piktogramme

STANDARD



OPTION



KERN SET-101

ET Save-cloud à Safe internet data storage

Software EasyTouch

Save-Cloud SET-101

Category

Brand: Accessories KERN

Category: Software

Product Group: Software

Packaging & shipping

Delivery: 24 h

Shipping method: Parcel service

Piktogramme

STANDARD



OPTION



KERN SET-151

ET BottleWeigh â Professional weighing of spirits / bottles, stock management bar statistics

Software EasyTouch

BottleWeigh SET-15

Category

Brand:	KERN
Category:	Software
Product Group:	Software

Packaging & shipping

Delivery:	24 h
Shipping method:	Parcel service

Piktogramme

STANDARD



OPTION



По вопросам продаж и поддержки обращайтесь:

Алматы (7273)495-231	Калининград (4012)72-03-81	Омск (3812)21-46-40	Сыктывкар (8212)25-95-17
Ангарск (3955)60-70-56	Калуга (4842)92-23-67	Орел (4862)44-53-42	Тамбов (4752)50-40-97
Архангельск (8182)63-90-72	Кемерово (3842)65-04-62	Оренбург (3532)37-68-04	Тверь (4822)63-31-35
Астрахань (8512)99-46-04	Киров (8332)68-02-04	Пенза (8412)22-31-16	Тольятти (8482)63-91-07
Барнаул (3852)73-04-60	Коломна (4966)23-41-49	Петрозаводск (8142)55-98-37	Томск (3822)98-41-53
Белгород (4722)40-23-64	Кострома (4942)77-07-48	Псков (8112)59-10-37	Тула (4872)33-79-87
Благовещенск (4162)22-76-07	Краснодар (861)203-40-90	Пермь (342)205-81-47	Тюмень (3452)66-21-18
Брянск (4832)59-03-52	Красноярск (391)204-63-61	Ростов-на-Дону (863)308-18-15	Ульяновск (8422)24-23-59
Владивосток (423)249-28-31	Курск (4712)77-13-04	Рязань (4912)46-61-64	Улан-Удэ (3012)59-97-51
Владикавказ (8672)28-90-48	Курган (3522)50-90-47	Самара (846)206-03-16	Уфа (347)229-48-12
Владимир (4922)49-43-18	Липецк (4742)52-20-81	Саранск (8342)22-96-24	Хабаровск (4212)92-98-04
Волгоград (844)278-03-48	Магнитогорск (3519)55-03-13	Санкт-Петербург (812)309-46-40	Чебоксары (8352)28-53-07
Вологда (8172)26-41-59	Москва (495)268-04-70	Саратов (845)249-38-78	Челябинск (351)202-03-61
Воронеж (473)204-51-73	Мурманск (8152)59-64-93	Севастополь (8692)22-31-93	Череповец (8202)49-02-64
Екатеринбург (343)384-55-89	Набережные Челны (8552)20-53-41	Симферополь (3652)67-13-56	Чита (3022)38-34-83
Иваново (4932)77-34-06	Нижний Новгород (831)429-08-12	Смоленск (4812)29-41-54	Якутск (4112)23-90-97
Ижевск (3412)26-03-58	Новокузнецк (3843)20-46-81	Сочи (862)225-72-31	Ярославль (4852)69-52-93
Иркутск (395)279-98-46	Ноябрьск (3496)41-32-12	Ставрополь (8652)20-65-13	
Казань (843)206-01-48	Новосибирск (383)227-86-73	Сургут (3462)77-98-35	
Россия +7(495)268-04-70	Киргизия +996(312)-96-26-47	Казахстан +7(7172)727-132	